

## Native Acrylamide Gels (35 ml)

	<b>3.5%</b>	<b>5%</b>	<b>6%</b>	<b>8%</b>
30/0.8% acrylamide	4.1 ml	5.8 ml	7 ml	9.3 ml
10X TBE	3.5 ml	3.5 ml	3.5 ml	3.5 ml
H <sub>2</sub> O	27.4 ml	25.7 ml	24.5 ml	22.2ml
Degas 1-2 min				

Add 0.25 ml 10% ammonium persulfate

Add 30 microliters TEMED

Pour Gel and wait 20-30 minutes before removing gel comb

Run at 180 volts in 1X TBE Buffer

Steve Hahn

Last modified Fri, Oct 16, 1998